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Introduktion

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Sequence designs

Normalized ensemble defect:	34.0 %	To Analysis
RNA_seq_strand	GGGACUACACGAAGGACGGGUCCGUGUCAACAUCAGUCUGAUAAGCUACACGUUGAGUAG AGUGUGAGCGUGUAGUCCCGUGAUUCCAGCGAAGGGACGGUGCGGAGAGGAGAGCUGGAA UCAC	<input checked="" type="checkbox"/>
Normalized ensemble defect:	34.9 %	To Analysis
RNA_seq_strand	GGGAUACACUGAAGGACGGGUCCGUGUCAACAUCAGUCUGAUAAGCUACACGUUGAGUAG AGUGUGAGCAGUGUAUCCCUAGGUGAAUGCGAAGGGACGGUGCGGAGAGGAGAGCAUUCA CCUA	<input checked="" type="checkbox"/>

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DNA template production:

1. Dissolve gBlock DNA to 100 nM and primers to 10 μ M in nuclease-free H₂O or 1X Tris acetate-EDTA (TAE) buffer (note 9).
2. Prepare PCR mix for a 50 μ l reaction: 10 μ l 5X Q5 Reaction buffer, 1 μ l 10 mM dNTP mix, 2.5 μ l of each 10 μ M forward and reverse primers, 2 μ l of 100 nM DNA template, 31.5 μ l nuclease-free H₂O.
3. Add 0.5 μ l of Q5 DNA polymerase, mix by pipetting and place the samples into the thermocycler, use a standard PCR procedure such as follows:
 - a. Initial denaturation at 98° C for 2 minutes
 - b. 30 cycles of:
 - i. Denaturation at 98°C for 10 seconds.
 - ii. Annealing at 69°C for 15 seconds (note 8).
 - iii. Polymerization at 72 °C for 20 seconds.
 - c. Final extension at 72°C for 2 minutes.
 - d. Storage at 4°C.
4. Purify using a standard PCR cleanup kit (Macherey-Nagel) following the manufacturer's instructions (note 10).

RNA production and purification:

5. DNA template production as stated in the previous section (3.1.2).
6. Prepare transcription by mixing DNA template, 1X transcription buffer 2, 1 mM. DTT, 2.5 mM each NTP, T7 RNA polymerase.
7. Incubate reaction at 37°C for 4 hours (note 13).
8. Stop reaction by adding 1 U/100 μ l of DNase I and incubating at 37°C for 15 mins.
9. Mix 1:1 with denaturing loading buffer and heat at 95°C for 5 mins.
10. Run at 15 W for 35 minutes.
11. Visualize gel bands using UV shadowing and cut the RNA bands out of the gel with a clean scalpel.
12. Elute overnight with nuclease-free H₂O.
13. Transfer the mix into the Freeze'n'Squeeze gel extraction spin columns to separate the liquid from the gel (note 14). Spin at 13000 x g for 3 minutes at room temperature.
14. Add 1/10 of the solution volume of 3M NaOAc to a final concentration of 0.3M NaOAc.

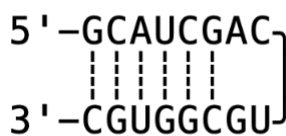
15. Add 3 times of the solution volume of 96% Ethanol, vortex and spin down.
16. Incubate sample on dry ice (-80°C) for 15-60 minutes or in the freezer (-20°C) for 1 hour to overnight.
17. Centrifuge at 14,000 rpm for 30 minutes. This should be done at 4°C if possible.
18. Carefully discard the supernatant and add 1.5 times the initial volume of 70% EtOH.
19. Centrifuge again at 14000rpm for 10 minutes.
20. Discard the supernatant and dry the pellet by leaving the tube in the fume-hood for a few minutes.
21. Re-dissolve the pellet in nuclease-free H₂O and quantify the concentration at the Nanodrop spectrophotometer using absorption at 260 nm. Store at -20°C.

Fluorescens-måling med RNA sensor og forskellige prøver:

22. Heat-anneal the RNA in the thermocycler by heating it up to 95 °C for 5 min, quickly cool it down by putting it in the freezer (-20°C) for 3 mins.
23. Add Assembly buffer to a final concentration of 1X and leave the samples at 37°C for 30 mins.
24. Prepare 100 nM samples for the fluorometer measurements.
25. Add 1 μM of DFHBI-1T and 1.7 μM YO3-biotin.
26. DFHBI-1T should be excited at 450 nm and its emission should be recorded at 503 nm and YO3-biotin should be excited at 580 nm and its emission should be recorded at 620 nm.
27. Relative FRET calculations should be done as in Jepsen *et al.* [10], by using the equation; $FRET = \frac{IDY(exD,emY)}{IDY(exD,emY)+IDY(exD,emD)}$, where $IDY(exD,emY)$ is the emission at YO3-biotin wavelength after DFHBI-1T excitation and $IDY(exD,emD)$ is the emission at DFHBI-1T wavelength after DFHBI-1T excitation. Both measurements should be performed with both fluorophores present.

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UACGCACGACUGGUAGCCCCUGGCACAGCCAGGUCCAGGAAAAAAGGA
UCAUGCGACGCAUGAGGCUACUUCCCUAUACGUGGUAUAGGGGUCGUG
CGC

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